Science with Passion

 **KNAUER**

*Azura*

**Organizer**

Supplement/Beiblatt

Document No. V6868

**Note:** Please read the corresponding technical documentation for handling and safety reasons.

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Version number: 2.0

Release date: 2022/12/28

The information in this document is subject to change without prior notice. For the latest version of the instructions, visit our website: [www.knauer.net/library](http://www.knauer.net/library).

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---

# Table of contents

|   |          |
|---|----------|
| <b>1. Product information .....</b>                             | <b>1</b> |
| <b>2. Scope of delivery .....</b>                               | <b>3</b> |
| <b>3. Attaching AZURA® Organizer onto AZURA Click rail.....</b> | <b>4</b> |
| 3.1 Attaching AZURA Organizer .....                             | 4        |
| 3.2 Taking off AZURA Organizer .....                            | 4        |
| <b>4. Mounting the clamps .....</b>                             | <b>5</b> |
| <b>5. Mounting AZURA pH flow cell .....</b>                     | <b>6</b> |
| <b>6. Repeat orders.....</b>                                    | <b>8</b> |

# 1. Product information

With the AZURA® Organizer, accessories such as columns, collection vessels, back pressure regulators or pH sensors can be grouped and fixed using a simple clamping mechanism.

## Legend:

- ① AZURA Organizer
- ② AZURA Click rail
- ③ AZURA device
- ④ Holes to mount the clamps
- ⑤ Holes to mount the AZURA pH flow cell
- ⑥ AZURA Click rail holder

Fig. 1 AZURA Organizer front side

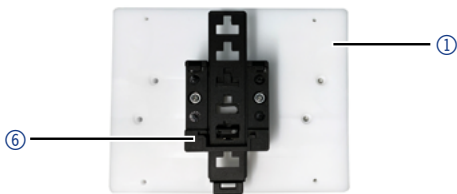
Fig. 2 AZURA Organizer back side

The AZURA Organizer can be flexibly positioned on the side surface of an AZURA device or system with L-footprint via the AZURA Click rail.

### Legend:

- ① AZURA Organizer
- ② 1 ml column
- ③ 5 ml column
- ④ AZURA pH flow cell
- ⑤ AZURA Click rail
- ⑥ 50 ml falcon tube
- ⑦ 15 ml falcon tube

**Fig. 3** AZURA Organizer with accessories

## 2. Scope of delivery

The delivery includes:

| Item               | Quantity | Usage  |
|--------------------|----------|--|
| AZURA Organizer    | 1        | to group and fixate clamps and accessoires, respectively   |
| 12 mm clamp        | 6        | for columns up to 13 mm outer diameter, e.g. 1 ml FPLC cartridges  |
| 16 mm clamp        | 4        | for 15 ml falcon tubes,<br>for columns up to 16 mm outer diameter, e.g. 5 ml FPLC cartridges or 120 ml SEC columns,<br>for a back pressure regulator |
| 25 mm clamp        | 2        | for 50 ml falcon tubes,<br>for columns up to 26 mm outer diameter  |
| Pan head screw T10 | 6        | to attach the clamps to the AZURA organizer  |

### Legend:

- ① AZURA Organizer
- ② Pan head screw T10 (6x)
- ③ 25 mm clamp (2x)
- ④ 16 mm clamp (4x)
- ⑤ 12 mm clamp (6x)

**Fig. 4** AZURA Organizer scope of delivery


## 3. Attaching AZURA® Organizer onto AZURA Click rail

### 3.1 Attaching AZURA Organizer

**Prerequisite:** The AZURA Click rail has been mounted.

#### Process

1. Hook the holder ① of the AZURA Organizer on the upper edge of the AZURA Click rail ②.
2. Push the AZURA Organizer against the lower edge of the AZURA Click rail.

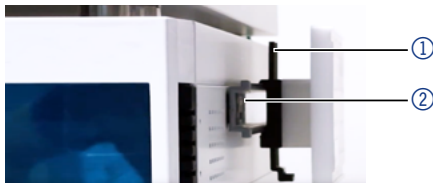**Fig. 5** Attaching AZURA Organizer on mounting rail

### 3.2 Taking off AZURA Organizer

#### Process

1. Pull the handle of the holder down.
2. Take off the AZURA Organizer by tilting it upwards.

**Video:** On the KNAUER YouTube channel you will find a short tutorial on how to attach and take off the AZURA Organizer from the mounting rail: [AZURA Click with AZURA Organizer](#).

## 4. Mounting the clamps

**Note:** When mounting the clamps:


- Use two clamps to mount longer columns.
- For SEC columns, KNAUER recommends two organizers.
- Use the screws supplied with the delivery.

**Tools:** ■ Torx screwdriver T10

- Pan head screws

### Process

1. Place the clamp ② on of the bore holes ③ of the AZURA Organizer.
2. Fasten the clamp with the pan head screw ① using the Torx screwdriver.
3. You can adjust the span of the clamp when tightening the screw.

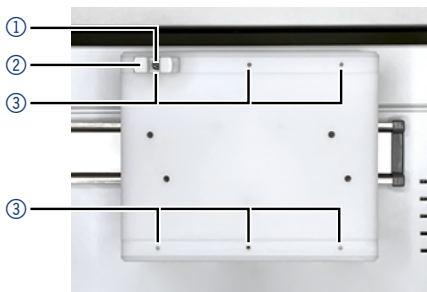**Fig. 6** Clamps on AZURA Organizer

## 5. Mounting AZURA pH flow cell


**Note:** In order to mount the AZURA pH flow cell onto the AZURA Organizer, the universal AZURA Click rail holder ① has to be unscrewed first.

- Tools:**
- Phillips screwdriver, PH1
  - Torx screwdriver, T10
  - 2 screws, size M3x06
  - 2 screws, size M3x10

### Process

1. Loosen the two screws ① on the back of the assembly with a Phillips screwdriver (PH1).
2. Remove the universal AZURA Click rail holder ②.

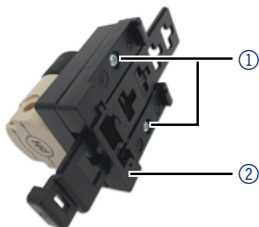**Fig. 7** Loosen two screws


3. Remove the AZURA Organizer ③ from the AZURA Click rail ④.

**Fig. 8** AZURA Organizer on AZURA Click rail (the holes for mounting the AZURA pH flow cell are each marked with a square and a circle)


## Process

- On the back side of the AZURA organizer **③** (where the AZURA Click rail holder **⑤** is attached), insert one of the screws **⑥** (size M3x10) into one of the holes (marked with a square/circle in the figure).

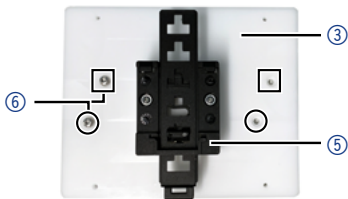**Fig. 9** Screws for fixing the AZURA pH flow cell on AZURA Organizer

- Hold the back of the AZURA pH flow cell **⑦** against the front of the AZURA Organizer so that the screw can be screwed into the corresponding threaded hole on the AZURA pH flow cell (marked with a square/circle in the picture). Use the T10 torx screwdriver. Turn the screw a few turns, but do not tighten it.

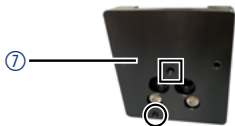**Fig. 10** Back of AZURA pH flow cell with corresponding threaded holes.

- Screw the the second screw into the other hole. Use the T10 torx screwdriver to tighten both screws slightly.

**Note:** It is sufficient to lightly tighten the screws to fix the AZURA pH flow cell to the AZURA Organizer.

## Process

- Attach the AZURA Organizer ③ with the mounted AZURA pH flow cell ⑦ to the AZURA Click rail ④.

**Fig. 11** AZURA Organizer with mounted AZURA pH flow cell on AZURA Click rail

**Note:** Alternatively, you can also attach the AZURA pH flow cell on the AZURA Organizer at the position of the other two holes.

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All rights reserved.
AZURA® is a registered trademark of KNAUER Wissenschaftliche Geräte GmbH.
Wissenschaftliche Geräte GmbH